



Immunofluorescence (IF) Protocol

Immunofluorescence (IF) microscopy is used to localize specific constituents in tissue sections or immobilized cells using fluorescent tags as labels. The antibody/antigen complex is labeled with any of a variety of fluorochromes emitting light from the near UV to the near IR. Usually FITC, a fluorescein derivative, is used for routine work and shows a yellow/green color. Red is often visualized using Texas Red, rhodamine, or Cy3. Blue occurs when AMCA or Cascade Blue is used. Fluorescent microscopes contain limit filters that are optimized for each fluorochrome. Confirm that your microscope is optimized for each dye you plan to use in your experiment. Black and white or color photography or enhanced imaging are often used to record data.

I. Reagents Required

Reagent	Preparation
Phosphate Buffered Saline (PBS) Wash Buffer (MB-008)	Use 10X PBS, pH 7.2 (0.2 M Potassium Phosphate, 1.5 M NaCl). Dilute appropriate volume to 1X with deionized water.
Formaldehyde Fixative	Dilute to 1% in PBS buffer. Use high quality formaldehyde.
Permeabilization Buffer	Prepare appropriate volume of 0.2% (v/v) Triton X-100 in PBS buffer.
Antibody Dilution Buffer	Prepare 100 mL of PBS Wash Buffer supplemented with 1 mL of normal serum of same species as the host used for the secondary antibody.
Fluorescein Secondary Antibody	Prepare dilution of fluorescein Conjugated secondary antibody (1:200) in Antibody Dilution Buffer. Use secondary antibody conjugate against the same species as the primary antibody.
Bis-Benzimide Solution	Prepare reagent at 0.5 mg/mL in PBS buffer.
Aqueous Mounting Media	N/A

II. Procedure

1. Grow cells on sterile 12 mm glass coverslips placed in 24-well culture plates. Remove culture medium.
2. Gently wash cells 3 times with ice cold PBS for 5 minutes per wash.
3. Fix cells by adding a volume of 1% formaldehyde in PBS equal to the original volume of culture medium. Incubate on ice for 5 minutes. Remove the fixative and wash as in Step 2.
4. Permeabilize cells for 20 minutes on ice with permeabilization buffer.
5. Prepare primary antibody appropriately diluted in antibody dilution buffer. React cells with primary antibody for 1 hour at room temperature. Wash as in Step 2.
6. Prepare dilution of fluorescein-conjugated, affinity-purified, secondary antibody. React cells for 1 hour at room temperature with reagent. Alternatively, use a biotin-conjugated secondary antibody followed by a wash as in Step 2.
7. Add fluorescein-conjugated streptavidin (S000-02) diluted 1:200 in PBS buffer. React for 30 minutes at room temperature. Wash as in Step 2.
8. Counterstain cells with bis-benzimide solution for 15 minutes at room temperature. Wash as in Step 2.
9. Add aqueous mounting agent. Affix coverslips to slides. Allow coverslips to dry in the dark before viewing.

References

- Ascoli and Maul. 1991. J. Cell Biol. 112:785-795.
- Antibody Techniques. V.S. Malik and E.P. Lillehoj, eds. Academic Press, 1994.

Rockland products are for research use only and are not intended for therapeutic or diagnostic applications. Please contact a technical service representative for more information. All properties listed are typical characteristics and are not specifications. All suggestions and data are offered in good faith but without guarantee as conditions and methods of use of our products are beyond our control. All claims must be made within 30 days following the date of delivery. Suggested uses of our products are not recommendations to use our products in violation of any patent or as a license under any patent of Rockland Immunochemicals, Inc. If you require a commercial license to use this material and do not have one, then return this material, unopened to: Rockland Immunochemicals, Inc., P.O. BOX 5199, Limerick, Pennsylvania 19468, USA.

© 2018 Rockland Immunochemicals, Inc. All rights reserved.