

# Immunohistochemistry (IHC) Protocol

## I. Reagents and Equipment Required

#### Reagent/Equipment

Phosphate Buffered Saline (PBS)(MB-008)

Xvlene

95% and 100% Ethanol

UltraPure Sterile Water (MB-009-1000)

Antibody Dilution Buffer: Prepare 100 mL of PBS, supplemented with 1 mL of normal serum of same species as host used for the secondary antibody.

30% Hydrogen Peroxide Solution (KHJ001)

Biotinylated Secondary Antibody, 1:500 (e.g. 611-106-122)

Streptavidin Peroxidase Conjugated, 1:500 (S000-03)

DAB Substrate (DAB-10) or TMB Membrane Peroxidase Substrate (TMBM-100) for stable brown or blue staining, respectively.

Polymount Mounting Media (KHH001)

#### II. Procedure for Frozen Sections

- 1. Snap-freeze fresh tissues in liquid nitrogen or isopentane pre-cooled in liquid nitrogen, embedded in OCT compound in cryomolds. Store frozen blocks at -80°C.
- Cut cryostat sections 4–8 mm thick and mount onto Superfrost™ Plus slides or gelatin-coated slides. Store slides at -80°C until needed.
- 3. Before staining, warm slides at room temperature for 30 minutes and fix in ice-cold acetone for 10 minutes. Air-dry for 30 minutes.
- 4. Wash in PBS.

### III. Procedure for Paraffin Sections

- 1. Deparaffinize sections in xylene 2 times for 5 minutes.
- 2. Hydrate with 100% ethanol 2 times for 3 minutes.
- 3. Hydrate with 95% ethanol for 1 minute.
- 4. Rinse in UltraPure sterile water.

## IV. Procedure for Immunoenzyme Staining

- 1. Follow procedure for pretreatment as required.
- 2. Rinse sections in PBS 2 times for 2 minutes.
- 3. Incubate sections in normal serum block with the same species as the secondary antibody (e.g., Normal Goat Serum (NGS) (B-304) if secondary antibody is goat host).
  - **Note:** Since this protocol uses avidin-biotin detection system, avidin/biotin block may be needed based on tissue type. If you do, the avidin/biotin blocking should be done after normal serum block and before primary antibody incubation.
- 4. Incubate sections in primary antibody at appropriate dilution in dilution buffer for 1 hour at room temperature or overnight.

  Note: Do not rinse sections between serum block and primary antibody incubation.
- 5. Rinse in PBS buffer for 3 times for 2 minutes.
- 6. Incubate sections in 1% hydrogen peroxidase in PBS for 10 minutes at room temperature.
- **7.** Rinse in PBS buffer 3 times for 2 minutes.
- 8. Incubate sections in biotinylated secondary antibody in PBS buffer for 30 minutes at room temperature.
- 9. Rinse in PBS buffer 3 times for 2 minutes.
- 10. Incubate sections in streptavidin peroxidase in PBS buffer for 30 minutes at room temperature.
- 11. Rinse in PBS buffer 3 times for 2 minutes.
- **12.** Incubate sections in peroxidase substrate solution.

- **13.** Rinse in PBS buffer 3 times for 2 minutes.
- 14. Rinse in UltraPure sterile water 3 times for 5 minutes.
- **15.** Dehydrate through 95% ethanol for 1 minute, 100% ethanol 2 times for 3 minutes.
- **16.** Clear in xylene 2 times for 5 minutes.
- 17. Coverslip with mounting medium (KHH001).

Rockland products are for research use only and are not intended for therapeutic or diagnostic applications. Please contact a technical service representative for more information. All properties listed are typical characteristics and are not specific ations. All suggestions and data are offered in good fairh but without guarantee as conditions and methods of use of our products are beyond our control. All relaterial and specifically into the product are product as a conditions and methods of use of our products are beyond our control. All relaterial and specific and pass following the death of delivery. Suggested uses of our products are not use our products are not used to the product are not used to the product are not used to the product are not product are not used to the product are not used t

© 2018 Rockland Immunochemicals, Inc. All rights reserved